



GOGGLE SANITIZER CABINET

S07817

Designed to meet the needs of science laboratories, classrooms, and factories. This heavy gauge CRC steel cabinet can sanitize up to 36 pairs of goggles in just 15 minutes using ultraviolet (UV) light. The goggle sanitizer cabinet is safe, easy to use, and can be placed on a desk, lab bench, or hung on the wall with the included mounting hardware.

FEATURES

- Made from heavy gauge CRC steel for durability with a white baked enamel finish for resistance to heat and corrosion.
- Internal metal racking system conveniently accommodates up to 36 goggles or safety-glasses with no stacking required.
- The front of the cabinet is equipped with a 15-minute timer, instructions, safety label, lock, and an on/off indicator light for safe and easy use.
- Tight piano style hinges, key lock, indicator light, and an automatic shut-off mechanism if the doors are opened during sterilization ensure everyone's safety.
- Cabinet operates with a long-life, high-intensity ultraviolet (UV) light bulb for reliable sanitizing and longevity.

OPERATION

Place eyewear in the cabinet, close and lock the doors. Set the timer for 15 minutes. The UV light will turn on and kill any germs present on the goggles, sterilizing them for the next user. The indicator light on the front of the cabinet will turn off when sterilization is complete and the UV bulb has been turned off.

SPECIFICATION

Input Voltage	120V AC \pm 10%, 50/60Hz.
Input Current	0.315A.
Wattage	25 Watts (Max).
Size	8.5 x 8.5 x 20 cm Deep
Weight	Approximately 54 lbs.
Cord Length	8 ft.
Inside Grid	3.5" x 3.5" x 8" Deep

In the United States:

For customer service, call 1-800-766-7000

To fax an order, use 1-800-926-1166

To order online: www.fishersci.com

www.fishersci.com/eisco