

HisTrap™ FF crude

Speed and convenience


HisTrap FF crude is a ready-to-use column for convenient purification of histidine-tagged proteins from unclarified lysates by immobilized metal ion affinity chromatography (IMAC).



- Reliable purification of histidine-tagged proteins directly from unclarified lysates – simply sonicate and run
- Reduced sample preparation time, which minimizes risk of protein degradation
- Compatible with a wide range of reducing agents, detergents, denaturants, and other additives
- Available in convenient, prepacked 1 ml and 5 ml HisTrap™ column formats
- Convenient HisTrap FF crude Kit offers three HisTrap FF crude 1 ml columns, premade buffers, connectors, and a syringe
- Simple operation with a syringe, pump, or chromatography system such as ÄKTA™ design


HisTrap FF crude

Enzymatic and mechanical cell lysis

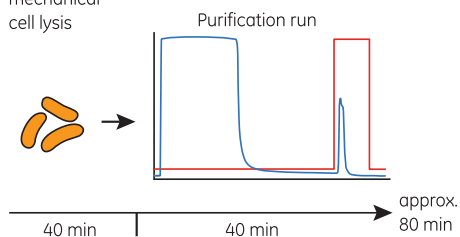

Conventional IMAC

Enzymatic and mechanical cell lysis

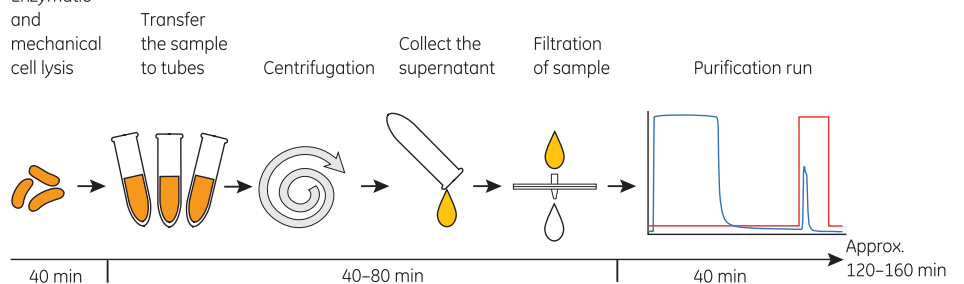

Schematic showing the time saved in purification of histidine-tagged proteins using HisTrap FF crude (only 80 min) compared to conventional IMAC (between 120 and 160 min).



HisTrap FF crude

Speed and convenience

HisTrap FF crude 1 ml and 5 ml columns permit rapid, yet reliable purifications from unclarified samples, with minimum sample preparation time and a reduced need for equipment.

Ordering information

Product	Quantity	Fisher Scientific Cat. No.
HisTrap FF crude	5 × 1 ml column	45000000
HisTrap FF crude	5 × 5 ml column	45000335
HisTrap FF crude Kit ¹	3 × 1 ml columns, premade buffers, connectors, syringe	45001428
His Buffer Kit ²	1 kit	45000007

¹ Contains 3 × 1 ml HisTrap FF crude columns, premade sodium phosphate buffer concentrate (2 × 50 ml phosphate buffer, pH 7.4), 2 M imidazole pH 7.4, connectors, syringe, and instructions.

² Contains premade sodium phosphate buffer concentrate (2 × 100 ml phosphate buffer, pH 7.4) and 2 M imidazole pH 7.4

To find out more about the complete Protein Purification product range visit www.gelifesciences.com/protein-purification

Your favorite Protein Purification products are now available from your local supplier.

GE, imagination at work, and GE monogram are trademarks of General Electric Company.

HisTrap is a trademark of GE Healthcare companies.

© 2011 General Electric Company—All rights reserved.

First published November 2009.



Distributor
GE Healthcare



In the United States:
For customer service, call 1-800-766-7000.
To fax an order, use 1-800-926-1166.
To order online: www.fishersci.com

In Canada:
For customer service, call 1-800-234-7437.
To fax an order, use 1-800-463-2996.
To order online: www.fishersci.ca