

NEULOG ACCELERATION LOGGER SENSOR GUIDE



NeuLog acceleration logger sensor NUL-227

The NeuLog acceleration sensor can be used for any science experiment or activity which utilizes acceleration measurements such as in the fields of Physics, Mechanics, Physiology, Exercise Science, etc.

The sensor comes pre-calibrated so you can start experimentation right out of the box using any of the following guides.

Just a few of the thousands of possible experimental subjects that can be done with the NUL-227 sensors are: gravitational force, impulse, exercise studies, human and animal acceleration, and many more


The acceleration sensor uses the following units of measure:

- Meters per second squared (m/s^2): The SI unit of acceleration or change in velocity over time.

Sensor orientation:

The NeuLog acceleration sensor has the ability to measure acceleration in three different axes (X, Y, and Z). The figure below demonstrates the axes.

Note: You can also easily determine which axis is which by using this easy method: First click on the sensor icon in the NeuLog software and select the axis you wish to determine. Then change the physical orientation of the acceleration sensor until the readout is roughly $\pm 9.8m/s^2$. When the sensor reads roughly $\pm 9.8m/s^2$, draw an imaginary line through the sensor towards the ground, either end of the imaginary line (sky and ground) will be the poles of the axis in relation to the sensor you are trying to determine.

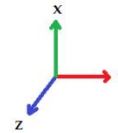

To view all axes simultaneously:

After collecting data:

1. Open the NeuLog software and connect the acceleration sensor using any of the following guides.
2. Once the sensor has been detected, click on "Offline Experiment"
3. Click on the acceleration sensor icon on the left.
4. Click on the "Range" button and select an axis (X, Y, or Z).
5. Click "Load Experiment" located on the top of the screen.
6. Select which experiment you would like to upload so that the button turns green, the default is "Newest" ("Newest" will select the most recent data collection)
7. Click on "Load Experiments" which will load the data from your chosen axis onto the graph.
8. Resume from step 3 to load the other axes on the same graph.

NEULOG ACCELERATION LOGGER SENSOR GUIDE



Quick start procedure:

PC or Mac Computer

Materials needed:

- NUL-227 Acceleration Sensor
- USB-200 USB Module
- A USB to mini USB cable (which comes with the USB-200)

Your acceleration sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via a USB to mini-USB cable. Please note that you cannot plug the acceleration sensor directly into the computer.

Resident PC software and browser based software can be downloaded for free at www.NeuLog.com/download as well as a full software user guide. **Note:** Make sure not to download and install both types of software, they will conflict on the computer.

Procedure:

1. Install the NeuLog software
2. Connect the USB-200 module to the PC or Mac
3. Connect the acceleration sensor to the USB-200 module (they directly plug together). Please note there is no calibration required for this sensor.
4. Open the NeuLog software.
5. Once a acceleration sensor logo appears on the left side of the screen the probe has been automatically identified and you can begin experimentation.
6. If the acceleration sensor is not automatically identified then click the "Search for sensors" icon to find the sensor.
7. Select the "On-line experiment" button; this will open a graph below.
8. Click on the "Module setup" button located on the acceleration sensor icon in the module window to change the sensor settings if need be.
9. Click on the experiment set up button to change the experiment settings if need be (experiment duration for example).
10. The acceleration sensor will give a live reading in the box to the left of the screen while plugged in.
11. To run an experiment and collect data click "Run experiment".
12. To end data collection early, click "Stop experiment".

Tablet, smart phone device

Materials needed:

- NUL-227 Acceleration Sensor
- WIFI-201 WIFI module
- BAT-200 Battery

Your acceleration sensor needs to be connected to a WIFI-201 module. The WIFI-201 module will create a closed NeuLog wifi network which will stream the NeuLog data to a device of your choosing. Once your device is wirelessly connected to the NeuLog network you can run experiments and collect data through a browser of your choosing.

Procedure:

1. Connect the acceleration sensor directly to the left side of a WIFI-201 module (no wires required).
2. Connect a BAT-200 module to the right side of the WIFI-201 module.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to micro USB charger (such as a typical cell phone charger). The WIFI-201 module will run for 15-75 minutes (depending on the sensor) without being plugged in.
4. For further WIFI-201 instructions or the WIFI-201 quick start guide please visit www.NeuLog.com/download.
5. The WIFI-201 indicator lights will flash; take no action until the LED to the far left turns blue.
6. Take your tablet or smart phone and go to the Wi-Fi settings and select the NeuLog network which matches the WIFI module ID found on the back of the WIFI-201 device. (NeuLogXXXX for example)
7. Give your device 1-2 minutes to connect to the WIFI-201 network.
8. Once the device is connected go to your browser and type in the website wifi201.com into the URL bar, then wait for 30-60 seconds.
9. You will see a "Control mode" icon in the browser, click on this icon.
10. The browser will then load a new screen and begin to auto detect the sensors, this can take a few minutes.
11. If the browser does not auto detect, select "Search for sensors"

NEULOG ACCELERATION LOGGER SENSOR GUIDE



12. Once the sensor is found you will see an icon on the left side of the screen for the acceleration sensor. The icon will display data in real time.
13. To set the acceleration sensor's settings click on "Module setup" button located on the acceleration sensor icon on the left side of your screen.
14. To change the experimental settings click the "Experiment setup" button (experiment duration for example).
15. To run an experiment and collect data click "Run experiment".

Operation with NeuLog viewer

Materials needed:

- NUL-227 Acceleration Sensor
- VIEW-101 Viewing Module
- BAT-200 Battery

Procedure:

1. Connect the acceleration sensor to the left side of the viewer.
2. Connect the BAT-200 module to the right side of the viewer.
3. The VIEW-101 will auto detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data click the run button (little green person).

Off-line experiments

(Off-line experiments are for when you do not have a sensor connected directly to a computer, tablet, smartphone, or NeuLog viewer).

Materials needed:

- NUL-227 Acceleration Sensor
- BAT-200 Battery

Materials needed to configure your offline experiment:

- USB-200 USB Module or WIFI-201 WIFI module or VIEW-101 Viewing Module
- A USB to mini USB cable (which comes with the USB-200)

Procedure:

1. Connect the acceleration sensor directly to the left side of a charged BAT-200 module.
2. When ready to collect data press the "Start/Stop" button on the acceleration sensor, a red light will turn on for the duration of the experiment.
3. To change the experiment settings first connect to your chosen device (PC, Mac, tablet, smart device) and select the "Module setup" button. (For more information on how to change the experiment settings view quick start procedure section for your chosen device in this document.)
4. After the experiment has concluded (when the "Start/Stop" button is pressed again or the data collection period ends) connect the sensor to your chosen device. (full instructions in each device's quick start procedure section)
5. Open NeuLog software.
6. Click the "Off-line experiment" button.
7. Click the "Load data from sensors" button.
8. Select which experimental data to upload. (5 experimental runs can be stored at one time on the NUL-227 acceleration sensor)

Restoring sensor's factory default settings:

Procedure:

1. Connect the USB-200 to a PC, Mac, tablet, or smart device.
2. Connect the acceleration sensor to a USB-200 module (they directly plug together). Please note there is no calibration required for this sensor.
3. Open the NeuLog software.
4. Click the "Tools" icon.
5. Click "Restore sensor's factory defaults" icon.
6. If prompted to clear the graph, click either the "Off-line experiment" or "On-line experiment" (whichever you are using).
7. Click "Clear experiment results".

NEULOG ACCELERATION LOGGER SENSOR GUIDE



8. Resume from step 4.

Included with sensor:

- NeuLog NUL-227 Acceleration Sensor Guide (this document).

Sensor specifications

Range and operation modes	$\pm 80 \text{ m/s}^2$
ADC resolution	10 bit
Resolution	0.15 m/s^2
Max sample rate (S/sec)	100 long duration 3000 short duration

Sensor features:

- Fully digital data.
- Rugged plastic ergonomic case.
- Push button switch for Start/Stop experiments in off line mode.
- Three internal accelerometers for X, Y, and Z axes.
- LED indicator of experiment status (blinks while collecting data).
- Pre-calibrated sensing equipment.
- **Note:** NeuLog products are intended for educational use.

Videos and experiment examples:

- Videos, literature and other probes can be found at www.NeuLog.com
- <http://www.NeuLog.com/Products/AccelerationloggerSensor.aspx>

Technical background:

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller

in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units ($^{\circ}\text{C}$, $^{\circ}\text{F}$, Lux, %, ppm, for example).

The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update.

The acceleration sensor houses three small thin silicon accelerometers, each accelerometer is oriented along a different axis (X, Y, or Z). When an accelerometer experiences an acceleration along its axis the silicon will flex to a very specific angle proportional to the acceleration it is experiencing. Each accelerometer is connected to a capacitor and as the silicon flexes the capacitance changes which in turn changes voltage flowing through the system. Acceleration and voltage have a linear relationship which allows the acceleration to be calculated easily using the voltage reading.

Maintenance and storage:

- Never submerge the NeuLog plastic body in any liquid.
- Do not allow liquid into the acceleration sensor's body.
- After use, gently wipe away any foreign material from the acceleration sensor.
- Store in a box at room temperature out of direct sunlight.

Warranty:

We promise to deliver our sensor free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your sensor will be repaired or replaced.

NEULOG ACCELERATION LOGGER SENSOR GUIDE



Thank you for using NeuLog!



Flexible, simple, fast, forward thinking.

W: www.neulog.com

E: info@neulog.com

A: 850 St Paul Street, Suite 15, Rochester, NY 14605

P: 1.866.553.8536